

committed to
brands.



**DURAN
WHEATON
KIMBLE**

Excellence in your hands

DURAN® **NEW** GL 25 AND GL 45 BROMOBUTYL RUBBER STOPPERS FOR EFFECTIVE SEALING

FEATURES

- Stopper shape, and material properties optimized to fit: 10 and 25 ml DURAN® GL 25 bottles, and DURAN® GL 45 bottles with sizes from 100 to 20,000 ml.
- **Low permeability** to air, gases and moisture.
- Manufactured from a **high purity bromobutyl rubber elastomer**.
- Elastic material for **effective sealing and reliable re-sealing** after needle piercing.
- Central needle piercing area indicated by mouled-in marks.
- Wide usable **temperature range: -50 to +121 °C**.
- Suitable for autoclaving / steam sterilizing at 121 °C for 20 minutes.
- Use with the **DURAN® GL 25 / GL 45 open topped / aperture screw cap** for extra security*.



BROMOBUTYL RUBBER – WHAT IS THAT?

Bromobutyl rubber is essentially impermeable to most gases, and provides a controlled environment inside the glass bottle for oxygen sensitive materials. Useful for maintaining anaerobic culture conditions. Bromobutyl rubber allows for multiple punctures providing easy access to the contents with a syringe.

HUBERLAB.

committed to science

DURAN® GL 25 AND GL 45 BROMOBUTYL RUBBER STOPPERS

TRUSTED QUALITY

- Conforms to **USP <381>** and **EP 3.2.9 Type I closure** for chemical and penetration tests.
- Material is virtually **free of organics** that might leach into the incubation medium.



APPLICATIONS

- Culture of **anaerobic microorganisms** (pressure resistant bottles recommended)
- Storage of **anaerobic culture media** (amber glass bottles recommended)
- Aerobic **methane oxidation (MOx)** research
- Storage of **oxygen** or **moisture sensitive** materials
- Applications requiring **gas purging** or **sample transfer** via a syringe

ORDERING INFORMATION

Description	Pack unit [pieces]	Cat. No.
NEW DURAN® Rubber Stopper for GL 25 laboratory bottles, straight plug, Grey Bromobutyl	10	29 206 13 01
NEW DURAN® Screw Cap GL 25 PBT red open topped with 15 mm central aperture	10	29 227 09 02
DURAN® Rubber Stopper for GL 45 laboratory bottles, straight plug, Grey Bromobutyl	10	29 206 28 03
DURAN® Screw Cap GL 45 PBT red open topped with 34 mm central aperture	10	29 227 10 07

Compatible products

Description	Pack unit	Cat. No.
DURAN® Laboratory Bottle, 10 ml, GL 25, clear	10	21 801 08 02
DURAN® Laboratory Bottle, 25 ml, GL 25, clear	10	21 801 14 04
DURAN® pressure plus+ Laboratory Bottle, GL 45, clear glass, 100 ml	10	21 810 24 06
DURAN® pressure plus+ Laboratory Bottle, GL 45, clear glass, 250 ml	10	10 922 34
DURAN® pressure plus+ Laboratory Bottle, GL 45, clear glass, 500 ml	10	10 922 35
DURAN® pressure plus+ Laboratory Bottle, GL 45, clear glass, 1000 ml	10	21 810 54 03
DURAN® pressure plus+ Laboratory Bottle, GL 45, amber glass, 100 ml	10	21 816 24 03
DURAN® pressure plus+ Laboratory Bottle, GL 45, amber glass, 250 ml	10	10 943 67
DURAN® pressure plus+ Laboratory Bottle, GL 45, amber glass, 500 ml	10	10 943 68
DURAN® pressure plus+ Laboratory Bottle, GL 45, amber glass, 1000 ml	10	21 816 54 09

* Standard DURAN® original GL 25 and GL 45 lab bottles are not pressure resistant. If build up in pressure is anticipated, it is recommended to use the stopper without the GL25 / GL45 red open top cap. The use of pressure resistant bottles such as DURAN® pressure plus+ should be considered.

001/3-2019



Contact.

HUBERLAB.

committed to science

OneStop. OneShop – HUBERLAB. bietet das Rundum-Sorglos-Paket für alle Bedürfnisse im Labor.

Profitieren Sie von unseren langjährigen Partnerschaften zu Top-Lieferanten.

Als Vollversorger für das Labor lässt unser umfassendes Sortiment keine Wünsche offen.

OneStop. OneShop – HUBERLAB. Nous offre le package sans souci pour tous les besoins de votre laboratoire.

Profitez de nos partenariats de longue date avec les meilleurs fournisseurs.

En tant que fournisseur complet pour le laboratoire, notre gamme complète ne laisse rien à désirer.

OneStop. OneShop – HUBERLAB. offers the all-included service package for all of your laboratory requirements.

Benefit from our long-standing partnerships with top-suppliers.

As a full-range supplier for the laboratory, our comprehensive range leaves nothing to be desired.



Wir beraten Sie gerne persönlich
Nous vous conseillons personnellement
We'll be happy to advise you personally



T +41 61 717 99 77



F +41 61 711 93 42



info.huberlab.ch



www@huberlab.ch



Let's chat
www.huberlab.ch



HUBERLAB. AG
Industriestrasse 123
4147 Aesch